

# Correlative Microscopy in Materials Science



Front cover: images from a workflow from X-ray microscopy (bulk inspection) to light microscopy (discovering an inclusion/particle) to electron microscopy (high-resolution image of the particle). Bottom right: complete additive manufactured part, non-destructively analysed with X-ray microscopy to check the severity of defects in general, then discovering a specific inclusion/particle with light microscopy (background image) and subsequent high-resolution imaging of the particle using electron microscopy for root cause determination (bottom left).

© 2017 John Wiley & Sons Ltd, The Atrium, Southern Gate,  
Chichester, West Sussex PO19 8SQ, UK  
Microscopy EKB Series Editor: Dr Julian Heath  
Spectroscopy and Separations EKB Series Editor: Nick Taylor

## CONTENTS

4 INTRODUCTION

6 HISTORY AND BACKGROUND

12 IN PRACTICE

24 PROBLEMS AND SOLUTIONS

29 WHAT'S NEXT?

### **About Essential Knowledge Briefings**

Essential Knowledge Briefings, published by John Wiley & Sons, comprise a series of short guides to the latest techniques, applications and equipment used in analytical science. Revised and updated annually, EKBs are an essential resource for scientists working in both academia and industry looking to update their understanding of key developments within each specialty. Free to download in a range of electronic formats, the EKB range is available at [www.essentialknowledgebriefings.com](http://www.essentialknowledgebriefings.com)

## INTRODUCTION

Correlative microscopy is not a single technique but a varied collection of techniques that share a common approach. By applying several different microscopy techniques to a single sample, scientists can study it at a much broader range of magnifications than is possible with a single technique. This allows them to conduct an initial low-magnification inspection of a sample to identify specific regions of interest (ROIs), which they can then zoom in on for a more detailed analysis, saving time and expense. There are other important advantages as well. Correlative microscopy allows scientists to study a greater diversity of samples, as some microscopy techniques work better with some materials than others, and to generate a much greater range of information about those samples at various different scales.

Combinations of light microscopy and electron microscopy are the most common form of correlative microscopy in use today. But researchers are increasingly designing correlative microscopy workflows that combine a greater number and variety of microscopy techniques, including X-ray microscopy, atomic force microscopy (AFM) and helium ion microscopy (HIM). These techniques are being combined to produce two-dimensional (2D) and three-dimensional (3D) images of a sample; by monitoring how samples change over time they are even producing four-dimensional (4D) images.

The demand for an ever-greater variety of correlative microscopy is being driven by the increasing range of materials that scientists want to study, and by the growing need to understand how the properties of those materials vary at different scales. While correlative microscopy was first applied to the study of

biological materials, scientists are now finding ways to apply it to a whole host of non-biological materials. These include geological materials, advanced alloys and nanomaterials such as graphene, as well as the advanced materials used in devices such as light-emitting diodes (LEDs), thin-film solar cells and batteries.

Microscope manufacturers are responding to this demand by developing ways to make correlative microscopy a more seamless and automated process. This includes developing software for combining the images, data and information generated by different microscopy techniques, advanced methods for identifying identical regions in images taken at different magnifications, and systems that physically combine different microscopy instruments (see Figure 1 on page 7).

This EKB offers an introduction to correlative microscopy. It begins by describing the different techniques that are commonly combined in correlative microscopy workflows, and outlines what specific benefits these different combinations provide. This is followed by a description of the combinations scientists tend to employ for studying different materials, with a specific focus on non-biological materials. There are also several case studies detailing specific examples of how scientists have applied correlative microscopy and what it has allowed them to discover.

The EKB also discusses the not-inconsiderable challenges of combining different microscopy techniques that work in different ways, require the sample to be treated in different ways and produce images at widely differing magnifications and resolutions. It finishes by highlighting some of the latest developments, including increasing automation and the introduction of super-resolution microscopy.

## HISTORY AND BACKGROUND


Correlative microscopy can be used to study an incredibly wide array of samples, from biological cells to multi-component technologies such as solar panels and lithium-ion batteries. Combining different microscopy techniques not only provides multi-scale insight (by allowing scientists to zoom in on a ROI) but also gives different types of complementary analytical information to maximize a study's usefulness. There are many cases where low magnification is required to survey large regions of a sample, for example to locate material defects, but high magnification is then required to obtain detailed structural information on those defects. Correlative microscopy allows researchers to meet both requirements.

This process is often assisted by the fact that light microscopy can produce color images of a sample, whereas high-magnification techniques such as electron microscopy can usually only produce black and white images. This means that ROIs that show up as specific colors, such as metal deposits, can be easily found in light microscopy images for more detailed study by electron microscopy, whereas they would be almost impossible to locate in electron microscopy images.

The ability to survey a material at low magnification in order to identify specific ROIs for study at higher magnifications offers several ancillary benefits. It allows scientists to use their microscopes more efficiently, both by locating ROIs faster at low magnification, thereby increasing analytical throughput, and by restricting the use of slower, more expensive, microscopy instruments to these identified ROIs. It also means that scientists aren't limited to the microscopes available in their own laboratory; samples can be transferred to another laboratory for study by different microscopes to reveal additional information.

Although scientists have long studied samples with different forms of microscopy, correlative microscopy as a discipline really began with the combination of light microscopy and electron microscopy to produce correlative light and electron microscopy (CLEM) (Figure 1). Commonly used in both life and material sciences, CLEM uses a light microscope to conduct a broad sweep of a sample, identifying ROIs that can then be studied in more detail with electron microscopy techniques such as scanning electron microscopy (SEM), focused ion beam (FIB)-SEM and transmission electron microscopy (TEM).

Correlative microscopy has since expanded to encompass many different microscopy techniques, covering a wide range of magnifications. Studies tend to begin with light microscopy, which has a maximum resolution of around 250nm. X-ray microscopy is the next step down: X-rays have shorter wavelengths than visible light and so X-ray microscopes can provide resolutions between those of light and

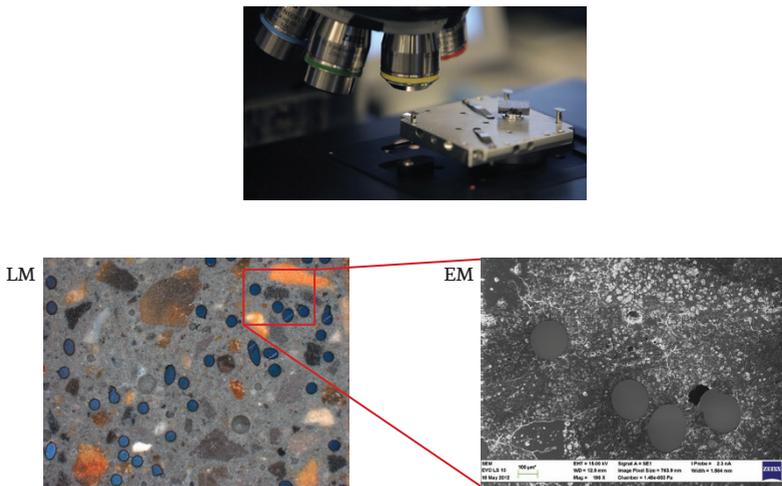

**Figure 1. Correlative microscopy combines microscopy techniques, for example light microscopy (LM) and electron microscopy (EM) to obtain maximal information about a sample. The top image shows a ZEISS flat samples holder for correlative microscopy**

electron microscopes, with the latest instruments able to achieve resolutions of up to 50nm. Furthermore, because X-rays penetrate most materials, performing X-ray microscopy on the same sample from various different angles allows researchers to build up 3D images of the sample. Known as X-ray computed tomography (CT), this process has a lower resolution than standard X-ray microscopy, able to achieve a maximum resolution of around 3.5 $\mu$ m. Furthermore, the achievable resolution tends to vary with the size of the sample.

TEM works by sending a beam of electrons straight through an ultra-thin sample section to a detector on the other side, with interactions between the electrons and the sample section generating the image. Because electrons have very short wavelengths, TEM can produce images at atomic-scale resolutions, able to show carbon atoms in diamond, for example.

SEM, on the other hand, scans an electron beam over a sample surface to generate secondary and backscattered electrons; these electrons can not only produce an image with a nanoscale resolution but can also reveal other information about the sample surface. Secondary electrons are released when the beam interacts with weakly-bonded electrons at the sample surface, and can reveal information about the topography (shape) and electrical properties of the sample. Backscattered electrons are electrons that have been reflected directly back from the sample. They tend to come from slightly deeper regions than secondary electrons and can reveal information about a sample's elemental make-up, because different atoms backscatter electrons to different extents.

Although SEM has a lower resolution than TEM, it does have several advantages. For instance, because it works by scanning an electron beam over a surface, it can analyze bulk samples, rather

than the thin sections required for TEM, and it also has a greater depth of view.

In addition to revealing information about a sample's topology and elemental make-up, the release of secondary electrons leaves behind positively charged 'holes' in the inner shells of atoms at the sample surface. As these holes fill with electrons from the outer shells, the difference in energy between the two shells is released in the form of an X-ray. Because the energies of these X-rays differ for different atoms, they can be used to determine the identity of the atom, in a technique known as energy-dispersive X-ray spectroscopy (EDS). Furthermore, the electron beam can also stimulate certain materials to emit light, in a process known as cathodoluminescence (CL). This can reveal information on the electronic structure of a sample and is particularly useful for determining the age of geological materials. All modern SEM instruments possess the detectors required for EDS and CL.


Many correlative techniques also make use of FIB-SEM, in which FIB and SEM are combined in a single system. FIB works in a similar way to SEM, but uses a beam of positively charged ions instead of negative electrons. This ion beam, which is usually produced by a liquid metal ion source such as gallium, can be used both to modify a sample and to analyze it. In an FIB-SEM system, the two instruments, which can be used simultaneously, are arranged so that the focal points of the electron and ion beam coincide.

At high ion doses, the FIB can dislodge atoms from the surface of a sample by sputtering, allowing it to modify or mill the sample. The FIB can also generate secondary electrons and ions, which can be used to produce images and provide information. For example, the secondary ions can be detected and identified by

secondary ion mass spectrometry (SIMS), revealing direct information about a sample's chemical composition. SIMS is more sensitive than EDS, able to detect chemical species at parts per million concentrations.

Using FIB-SEM in correlative studies allows researchers to look beneath a sample's surface. Once an ROI has been identified, the FIB can carve away successive layers of a sample while the SEM produces images of each newly uncovered surface. These images can then be combined to produce a 3D image of the ROI, in a process known as FIB-SEM tomography (Figure 2). FIB can also produce the thin sample sections required for TEM.

In some specialized cases, light microscopy can also be combined with HIM, which images samples with a beam of helium ions. The lighter helium ions do not penetrate as deeply into samples as the liquid metal ions used in FIB-SEM, producing secondary electrons and ions from the top 25nm of the sample surface.

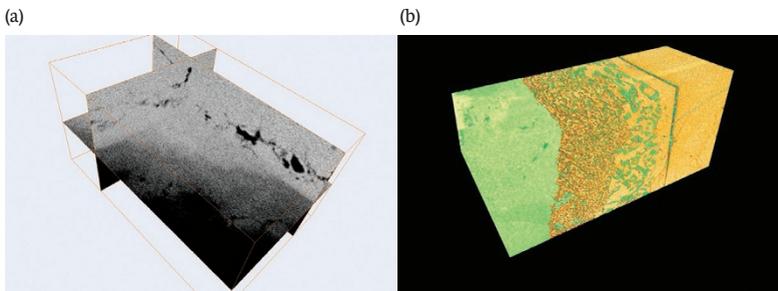

**Figure 2. (a) Volume rendering (yellow line) of X-ray microscopy dataset (approximately 0.5mm field of view) showing transition zone defects and coarse diffusion pathways in this thermoelectric system. 200nm voxel acquisition, showing three phases (black -  $\text{Mg}_2\text{Sn}$ , medium gray - transition zone, light gray -  $\text{Mg}_2\text{Si}$ ). (b) Volume rendering of ZEISS FIB-SEM acquisition. Silicon-rich phase is depicted in orange, and the stannum-rich phase in green, with clear interdiffusion in the transition zone approximately 25–30 $\mu\text{m}$  wide in the center**

Able to achieve a depth of field resolution of just 0.5nm – five to ten times greater than can be achieved with FIB-SEM – HIM has proved useful for studying very thin layers of graphene, as well as the thin films used in the electronics industry. It is also indispensable for studying samples containing gallium, which can be contaminated by a FIB made up of gallium ions.

‘Low dimensional’ materials such as graphene are also sometimes studied using Raman microscopy, which detects the scattering of light by molecular vibrations. For example, Raman imaging has been used to study strain and disorder in graphene. Because these molecular vibrations and associated light scattering are characteristic for specific molecules, Raman microscopy can also be used to determine the chemical composition of samples.

Correlative microscopy has already proved invaluable in two dimensions. With ROIs often lurking below the sample surface, often a few tens of micrometers below, many researchers are now also seeking 3D insights when studying networks of structures such as pore channels in rocks, fibers in composites or cracks in alloys. This so-called correlative tomography brings together multiple microscopy techniques able to study a 3D volume of interest, particularly X-ray microscopy and FIB-SEM tomography. Like correlative microscopy, correlative tomography provides access to a far larger range of scales for the same region than is possible with a single instrument, allowing scientists to investigate how material properties differ at different scales.

Running 3D experiments over time (time-lapse) adds a fourth dimension. Such 4D experiments are particularly valuable for monitoring changes in materials, such as following how cracks spread or spotting the moment at which metal fatigue occurs (see Case Study 2).

## IN PRACTICE

Correlative microscopy has already been utilized for a wide range of scientific disciplines, from biology, zoology and palaeontology, to materials and electronics. For example, a single team at the University of Manchester in the UK, led by Professor Philip Withers, has used correlative microscopy to study chrysalises in real time, structural changes in rats' arteries, metal fatigue and cracking, and woven composite materials such as wind turbine blades.

One of the great advantages of correlative microscopy is that, because it encompasses a range of different techniques, scientists can often find just the right combination for their studies. This is especially useful when studying non-biological materials that can differ widely in composition, structure and physical properties.


CLEM has proved invaluable for studying optoelectronic devices comprising different optical materials and photonic structures. For example, LEDs can be analyzed using light microscopy and SEM. Light microscopy provides a quick overview of morphology and optical properties such as color or reflectivity; the backscattered electrons generated by SEM can then be used to visualize contrast in cross sections, while EDS can determine elemental composition.

The same techniques are also being used to study thin-film solar cells, which are designed to provide a low-cost alternative to silicon wafer-based cells. The structure of the interfaces between different layers in these cells can be studied at an atomic scale with scanning TEM (STEM) while EDS can provide information on elemental distribution. In addition, SEM can provide insights

into the structural, compositional and electrical properties of the thin films, while backscattered electrons can give crystallographic information.

In geology, CLEM, together with CL, can be used to unearth a rock's age and deformation history. Meanwhile, 4D correlative microscopy can reveal how rocks 'corrode' over time.

A combination of X-ray microscopy and FIB-SEM has proved useful for studying lithium-ion batteries, as part of efforts to boost their performance and lifetime. In a study at Aalen University in Germany, X-ray microscopy revealed bright spots below the surface of a cathode from a commercial battery (Figure 3a). During subsequent FIB-SEM analysis, the researchers collected secondary electron and backscattered images that again showed

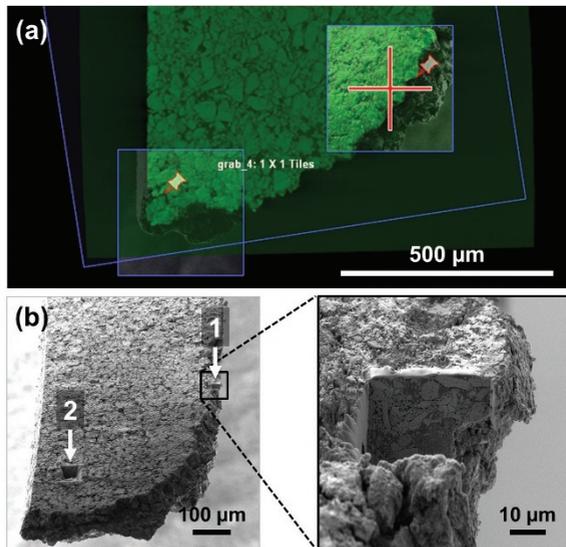

**Figure 3.** (a) Overlay of SEM (gray) and X-ray microscopy (green) images. After successful correlation, the X-ray microscopy data is used as a reference for FIB-SEM navigation. (b) SEM overview of the sample. Sites of interest are highlighted by the white arrows. The inset shows a detail of site 1

the particles (Figure 3b). EDS revealed these particles to be rich in lanthanum, rather than the expected manganese.

Other teams have used correlative microscopy to study the structure of batteries, for example how they contract and expand during charging. Over many charging cycles, batteries become more damaged, so understanding this process could help to develop longer-lasting products.

## Case study 1. Additive manufacturing

Additive manufacturing, also known as 3D printing, fabricates items layer-by-layer according to a computer-based design. It is perfect for producing engineering components, allowing parts with complex geometries and integrated functional properties, such as curved cooling channels, to be constructed from metal, alloy or ceramic powders.

As well as a raft of benefits, additive manufacturing also presents significant challenges, including the fact that 3D printed structures can be prone to forming microstructural defects such as small holes and weak areas. Correlative microscopy can help to root out these structural faults and understand why they are occurring.

At the Materials Research Institute in Aalen University, Germany, Tim Schubert, Dr Timo Bernthaler and Professor Gerhard Schneider have used correlative microscopy to find voids and inclusions in aluminum gear parts made using an additive manufacturing technique called laser powder bed fusion (see the images on the front cover of this *EKB*). For this, a laser melts a layer of metal powder deposited in the shape of a 2D slice through a 3D computer-based design. Once this layer has solidified, another layer of powder, representing the next 2D slice, is added on top, and the process repeated until the part is complete.

The team first used X-ray microscopy to conduct high-resolution imaging of the whole gear part, which allowed them to identify voids and inclusions. Next, they studied a micro-section of the gear with both light microscopy and SEM, using a purpose-built sample holder to ensure they were zooming in on the correct regions of the gear.

This meant they could study the inclusions identified by light microscopy with EDS, which revealed unexpectedly high levels of a titanium impurity. This titanium could be derived from impurities in the powder

raw material, suggests Bernthaler, or even from something as simple as insufficient cleaning.

‘More digitalization (software tools) and automation of microscopy have pushed correlative technology in recent years,’ says Bernthaler. By using ‘clever sample holders’ and software, workflows have greatly cut the time that such investigations take, he adds.

**Weissmayer L, Schubert T, Bernthaler T, et al. Applications of microscopy in additive manufacturing. *Optik & Photonik* 2015;4:44–6. (<http://dx.doi.org/10.1002/opph.201500015>)**

## Case study 2. Corrosion in magnesium alloys


Magnesium alloys are lightweight, high-strength alternatives to aluminum-based products for use in certain automotive and aerospace applications. But these alloys are sensitive to impurities, which can promote the spread of any damage, and are also prone to corrosion in humid or wet environments. With his team at Manchester University in the UK, Philip Withers, Regius Professor of Materials, has investigated how corrosion damage occurs in these magnesium alloys over time, at scales ranging from millimeters to nanometers and both at and below the surface.

To do this, Withers and his team first studied the alloy with X-ray microscopy while it was immersed in a corrosive salt solution, conducting a single analysis every two hours. In this way, Withers and his team produced a series of 3D images of the alloy corroding over time. They then studied the corroded alloy with SEM and combined this image with the 3D images produced by X-ray microscopy in order to identify a highly corroded ROI within the bulk of the sample.

After identifying such a region, they cut it out of the alloy with FIB to produce a microscopic sub-sample. This required relating the detailed surface image generated by SEM with the 3D X-ray images in order to determine where to remove material from the surface in order to reach the ROI. Next, they studied the sub-sample with X-ray microscopy, but using an X-ray microscope that could produce 3D images at a higher level of resolution. To increase the magnification further, they also produced a 3D image with FIB-SEM tomography. This involved removing multiple layers of the sub-sample with FIB and then imaging each revealed surface with SEM, before combining all the images together to produce the 3D image.

With X-ray microscopy, Withers and his team were able to see pitting across the magnesium alloy's surface within several hours of immersion in the salt solution. After 14 hours, the corrosion had spread below the surface, as revealed by FIB-SEM. By understanding how corrosion spreads, the researchers hope to find a way to prevent it.

Withers and his colleagues regularly use correlative microscopy to study cracks in engineering materials, which relate to defects, inclusions and other features. The key is to study the systems in 4D, says Withers. 'If you were to study lots of samples, including fatigued ones, it wouldn't necessarily be the case that you would get the same kind of crack growth,' explains Withers. 'But if you can study the same sample, we can actually watch how the cracks nucleate, how they connect up, and eventually lead to failure.'

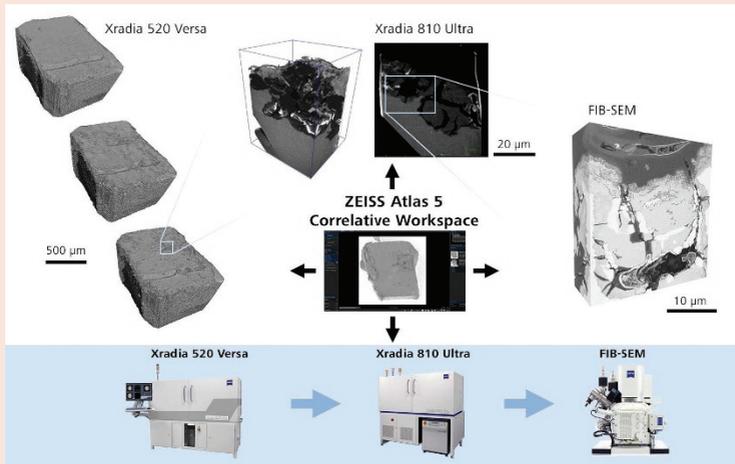

**Experimental overview.** This multiscale correlation study was designed to better understand corrosion damage in a Mg AZ318 alloy sample. The Atlas 5 correlative workspace was used to link and navigate between imaging performed on Xradia 520 Versa, Xradia 810 Ultra and the FIB-SEM microscopes

*In future, he predicts that further developments in automating the correlation between different instruments and imaging modalities will bring a ‘rapid expansion’ of multiscale correlative microscopy.*

**Harris, W. Multi-scale correlative study of corrosion evolution in a magnesium alloy. ZEISS Application Note, December 2015.**

### Case study 3. Studying structures in 3D over time

*‘We feel like kids in a candy store. We have all of these new tools. And we are really able to get inside problems that we haven’t been able to solve to date,’ says Nikhilesh Chawla, Fulton Professor of Materials Science and Engineering at Arizona State University in the USA.*

*Chawla is head of the new Center for 4D Materials Science (4DMS), funded by a collaboration between Arizona State University, ZEISS and the US Office of Naval Research. The center will use correlative microscopy to study structures in 3D over time.*

*‘The idea is to really push the state of the art in 4D material science,’ says Chawla. ‘We are essentially hoping to use these capabilities to solve problems in a variety of different areas, ranging from material science to civil engineering, mechanical engineering, electrical engineering and other physical sciences.’*

*Much of Chawla’s correlative microscopy brings together X-ray microscopy and electron microscopy, including FIB-SEM. His team uses ZEISS software called Atlas 5 to integrate data from the different techniques. ‘Atlas 5 really is revolutionary because in previous analyses finding the same region in the CT and in the FIB was like finding a needle in a haystack,’ he says.*

*For example, his team has applied non-destructive X-ray microscopy and FIB-SEM to the study of corrosion in aluminum alloys.<sup>1</sup> Using FIB to expose cross-sections of an alloy sample, perpendicular to the surface, while simultaneously using SEM to image the cross-section face revealed magnesium-rich precipitates inside the structure.*

*‘We were able to use compositional information to determine that the magnesium precipitates are actually dissolving preferentially in the corrosive*

fluid,' says Chawla. 'Other kinds of precipitates or particles based on iron, for example, were not really touched so that told us that the magnesium-rich precipitates are really the ones that are at risk in those kinds of alloys.'

Illustrating the diversity of his work, Chawla's team is also using correlative microscopy to study beetles called weevils,<sup>2</sup> which have very long and flexible 'beaks' for boring into plants. Using X-ray CT microscopy and FIB, the researchers have revealed that the beak structure comprises alternating hard and soft layers. 'The soft tissue is a kind of composite. It's unique. We can use the intelligent structures in the weevils to design new, man-made synthetic composites,' says Chawla.

Next, the researchers will incorporate a technique called nano-indentation, which employs a sharp diamond probe to test mechanical properties, into their workflow. 'That is the next avenue, to couple the structural information from CT or FIB with mechanical property information in site. That's something we can do inside the ion beam,' he explains.

1. Singh S, Loza J, Merkle A, et al. Three dimensional microstructural characterization of nanoscale precipitates in AA7075-T651 by focused ion beam (FIB) tomography. *Mater Charact* 2016;118:102-11. (<http://dx.doi.org/10.1016/j.matchar.2016.05.009>)
2. Andrew Jansen M, Singh, S, Chawla N, et al. A multilayer micro-mechanical model of the cuticle of *Curculio longinasus* Chittenden, 1927 (Coleoptera: Curculionidae). *J Struct Biol* 2016;195:139-58. (<http://dx.doi.org/10.1016/j.jsb.2016.05.007>)

## Case study 4. Polyurethane liners in splashdown airbags

One day, NASA's Orion crew module will carry astronauts beyond the moon, piggybacking on a new space launch system rocket. Orion's first uncrewed exploration flight took place in 2014, with the module safely splashing down in the Pacific Ocean. Unfortunately, however, the test revealed a failure in Orion's crew module uprighting system.

This system consists of five airbags that inflate upon splash-down, keeping the module upright until it can be rescued by the navy. Lab tests using correlative microscopy later revealed a problem with the polyurethane liner in the airbags, says James Martinez, engineering technologist at NASA's Lyndon B Johnson Space Center in Houston, Texas.

These tests suggested that the polyurethane layers were fusing or binding to each other under high pressure, causing them to tear. 'We are changing to another material that has a slip agent inside so that the two surfaces don't bond to themselves and can slip and slide against each other when the bag is deployed,' says Martinez.

When they studied this new layer with optical microscopy, however, Martinez and his team spotted inclusions in the polyurethane thin film. Areas of birefringence near the inclusions pointed to strain.

Using EDS, the team discovered that the slip agent in the material they had purchased is a crushed wafer-like substance called diatomaceous earth. Concerned that this might cause abrasion when the polyurethane layers rub against each other, the team used SEM at a low voltage (to prevent the material from degrading during imaging) to look for surface scratches or particulates. They then used SEM at a high voltage to study both the inclusions and the slip agent.

*‘It’s the best example of correlative microscopy that we have had in some time. The nature of this material really lends itself well to imaging with different modes and techniques,’ says Martinez. NASA is now carrying out a further series of safety tests, including on the new liner.*


## PROBLEMS AND SOLUTIONS

An effective workflow is essential for getting the most out of correlative microscopy. Such a workflow defines in what order a sample will be studied by the different microscopy instruments, how ROIs on the sample will be selected and identified in different instruments, and how the resultant data will be collected, stored and processed.

Usually, the level of magnification will steadily increase as the workflow progresses, in order to study a sample in progressively greater detail. But that's not always the sole reason for conducting correlative microscopy, and these other reasons may need to be reflected in the workflow.

For example, researchers may want to take advantage of the fact that different microscopy techniques can generate different information about a sample, which can be overlaid on the visual images. This is particularly the case with FIB-SEM, which can generate information about the elemental composition of a sample through both EDS and SIMS.

Alternatively, different microscopy techniques may be required to study the different materials in a sample, as certain materials may not respond well to specific techniques. For example, SEM works best with conducting substances such as metals, and both it and EDS can struggle with materials containing light elements such as carbon and lithium. This can clearly create problems when studying lithium-ion batteries, which employ lithium-based materials as the cathode and electrolyte, and graphite as the anode. One solution is to combine SEM with Raman microscopy (Figure 4), which can't analyze metals but can identify light elements and organic compounds.

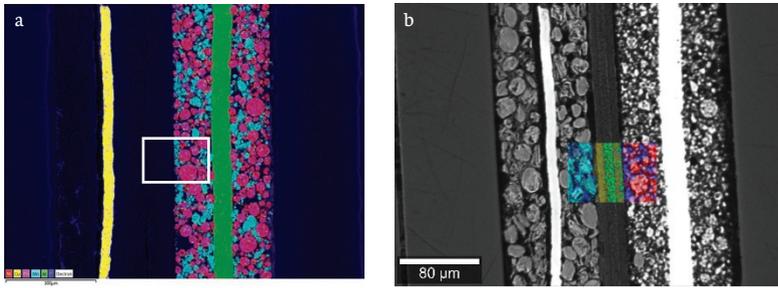

**Figure 4. (a) SEM image with EDS mapping. (b) Light microscopy and Raman overlay of exactly the same region**

Users also increasingly want to swap between 2D and 3D images produced by different techniques. Often they will do this merely to help identify a ROI for more detailed analysis, such as conducting X-ray microscopy to produce a 3D image that will reveal interesting features for closer study. But it can also involve conducting various rounds of 2D and 3D analysis with different techniques. Indeed, the development of correlative microscopy over the past few years has generally seen a progression from purely 2D workflows to those that incorporate more and more 3D techniques. Formerly the preserve of large universities, 3D correlative microscopy is now being adopted by smaller universities and industry.

One example is provided by the work of Professor Philip Withers and his team at Manchester University, which is detailed in Case Study 2. Another example is provided by a recent study conducted by Dr Paul Shearing and his team at University College London in the UK, who used correlative microscopy in 2D and 3D to study a commercial lithium-ion battery. This involved using X-ray microscopy to study the battery in 3D at a range of resolutions, from 22 $\mu\text{m}$  to 130nm, with two different microscopes, and FIB-SEM to reveal and study a sub-surface ROI in 2D by both SEM and EDS. This approach allowed Shearing and his team to study

the battery at a range of different scales, from the whole battery to the specific ROI, revealing information on everything from the large-scale structure of the battery to the nanoscale porosity of the electrodes.

This raises a critical factor that needs to be borne in mind when designing a workflow for correlative microscopy. Such workflows will often involve a combination of non-destructive techniques, such as X-ray microscopy, and destructive techniques, such as FIB-SEM, and clearly the position of any destructive technique in the workflow is critical. In particular, FIB-SEM tomography, in which successive layers are removed from a ROI to build up a 3D image, obviously needs to be conducted toward the end of a workflow, as it destroys the ROI. But any treatment or modification to a sample required for a specific microscopy technique may also influence the sequence of techniques.

Now, this is more of an issue when studying biological samples, which are often analyzed by light microscopy with minimal treatment but need to be encased in resin or flash-frozen prior to analysis by electron microscopy. While such treatments aren't required for non-biological samples, non-conductive materials may need to be covered in a thin layer of a metal such as platinum to improve the contrast in electron microscopy, which would influence the workflow. Such layers can also be applied to protect ROIs when milling them with FIB.

Once an appropriate workflow has been designed, users then need an effective way to collate and combine the data produced by the different microscopy techniques. Several software tools, such as Atlas 5 from ZEISS and MAPS from FEI, will automatically collect the data from a range of different techniques; for Atlas 5, this can

include various forms of light microscopy, X-ray microscopy and electron microscopy. They can then produce composite images of a sample by overlaying images produced by the different techniques at different magnifications, as well as overlaying chemical information produced by techniques such as EDS and SIMS.

For the software tools to be able to overlay images in this way, however, users need to match identical regions of the sample in the images produced by the different techniques. This can be done by marking the region in some way, such as by simply circling it with a pen, but this is not always desirable or possible. Alternatively, it can be done by identifying the same characteristic features in different images, but such features can look very different at different scales. Transferring the sample between different instruments, such that it is studied at slightly different orientations, adds further difficulty, as does any treatment or modification applied to the sample.

One way to make this process easier is to mount the sample on a holder that can be transferred between different microscopy instruments, thereby ensuring that the sample is always studied in the same orientation. Fiducial markers can also be added to the holder, providing fixed reference points that act as landmarks for locating the same regions and features in different images.

This is what ZEISS has done with its 'Shuttle & Find' technique, which employs a holder with three L-shaped fiducial markers that can be used with many of the light and electron microscopes produced by ZEISS. By measuring the distance from these markers to features on the sample, the Shuttle & Find software module can construct a detailed co-ordinate system for accurately and automatically locating the same region of the sample in different images.

This allows users to tag ROIs in images produced by a light microscope and then easily find those ROIs in images of the same sample produced by an electron microscope.

Some manufacturers have combined light microscopy and electron microscopy in single systems, removing the need to transfer samples between individual instruments and making it easier to match images of the same region produced by the two techniques. The downside to this approach is that not only does it restrict correlative microscopy to just these two techniques, but physically combining light and electron microscopy in a single system tends to compromise their performance. It also means that the two techniques can't be used to study different samples at the same time, reducing analytical throughput.

That just leaves the challenge of storing, processing and sharing all the data that can now be generated by correlative microscopy. This data is often substantial, especially if it includes 3D data sets, and thus requires lots of computer memory and processing power. Storing and processing this data in the cloud would appear to be an obvious solution, especially as it would also offer an easy way to share the data among different researchers. The problem with this approach is the great length of time it takes to upload all the data into the cloud. This data challenge is shared by many other areas of science, however, and solving it will require more than just advances in correlative microscopy.

## WHAT'S NEXT

Increased automation would help to resolve many of the current difficulties with correlative microscopy. Ideally, users would be able to study a sample on two or more instruments and then simply press a button to combine and correlate all the images, allowing the same region to be investigated seamlessly at a broad range of different scales.

Modern software tools such as Atlas 5, together with practical techniques such as Shuttle & Find, are going some way towards achieving this aim. Already, they are making it easier for researchers to analyze materials in two, three and even four dimensions and various different scales, and then combine the different sources of information. Nevertheless, other tools will be required to fully achieve this aim, including an automated method for revealing ROIs located within the bulk of a sample.

Researchers such as those profiled in the four case studies now regularly utilize these tools and techniques in their work. This is leading to the first practical applications of correlative microscopy, beyond purely academic research. For example, it is now being used for *in-situ* testing. This involves using a technique like X-ray microscopy to map the pores and voids in a material such as a metal alloy and then SEM to monitor the development of these pores and voids as the material is stressed until it breaks.

So far, these tools and techniques mainly work with light, X-ray and electron microscopes, but similar automated approaches are required for other microscopy techniques. Especially as researchers are starting to expand correlative microscopy to encompass a broader range of techniques.

One of these is AFM. Rather than produce a magnified optical image of a sample, AFM uses a flexible cantilever to measure the forces between the tip of the cantilever and the surface of a sample, as the tip is scanned across. Because these forces vary depending on the distance between the tip and the sample, they can be used to measure variations in the height of the surface, allowing a map of surface topology to be built up. Using the same basic approach, AFM can also probe a wide array of other material properties, including adhesion, elasticity, conductivity and temperature.

AFM is often combined with light microscopy, in order to compare the topological map with a surface image and also to direct the AFM analysis. As with light and electron microscopy, the problem is matching up images produced at widely different magnifications, as AFM is able to resolve features below 1nm size. The recent development of several super-resolution microscopy techniques, which were collectively awarded the Nobel prize for chemistry in 2014, offers a novel solution to this problem. These techniques offer ways to bypass the diffraction limit that normally limits the resolution of light microscopy to 250nm, bringing their resolution closer to that of AFM.

Up to now, the combination of AFM and super-resolution microscopy has mainly been used to study biological material, because super-resolution microscopy makes use of the fluorescent probes commonly used to illuminate cellular components in conventional fluorescence microscopy. Recently, however, a team of scientists led by Dr Cristina Flors at the Madrid Institute for Advanced Studies in Nanoscience in Spain used a combination of AFM and a localization-based super-resolution microscopy technique to study hybrid nanomaterials.

These hybrid nanomaterials comprised amyloid-like fibrils covered in the fluorescent nanocrystals known as quantum dots; such combinations of protein fibrils and nanoparticles are being developed for various applications, including as novel electrical switches and biosensors. The presence of fluorescent quantum dots on the hybrid nanomaterial made it ideal for study by super-resolution microscopy. Using the two techniques, Flors and her team were able to obtain useful information about both the structure of the amyloid-like fibrils and the distribution of quantum dots over the fibrils.

Super-resolution microscopy is also being combined with many other microscopy techniques, including light microscopy techniques such as fluorescence microscopy and bright-field microscopy, as an alternative to electron microscopy. Thus scientists can identify interesting features on a sample with light microscopy and then zoom in for a closer look with super-resolution microscopy. It is also being combined with electron microscopy, as both techniques have comparable resolutions.


As with AFM and super-resolution microscopy, these combinations have so far mainly been applied to the study of biological samples, but there have been a few exceptions. For example, a combination of SEM and localization-based super-resolution microscopy techniques have been used to study the structure of gold nanowires covered in fluorescent probes, as well as the activity of gold nanorods and gold nanoplates able to catalyze a reaction that produces a fluorescent product. For the gold nanorods, this revealed that their catalytic activity is higher at each end than in the middle.

Similar advanced nanomaterials are also being studied by other forms of correlative microscopy. These include graphene, a

single-atom-thick layer of carbon dubbed a ‘wonder material’ for its impressive properties, including great strength and extremely high electrical conductivity. However, graphene is often produced as flakes comprising several different layers, and the exact thickness of these flakes can have a major influence on their properties.

Scientists from ZEISS and DME Nanotechnologie (now Semi-lab) have shown how Shuttle & Find can be used to locate individual graphene flakes with a light microscope for more detailed study by SEM and AFM. This allowed the scientists to measure the precise thickness of different graphene flakes, and also revealed that the electronic and optical properties of graphene differ depending on the substrate it is placed on (Figure 5).

As graphene comprises atom-thick layers, it is an example of a 2D crystal. Another example is transition metal dichalogenides (TMDCs), a family of 2D crystals made up of a layer of transition metal atoms sandwiched between two layers of chalcogen atoms,

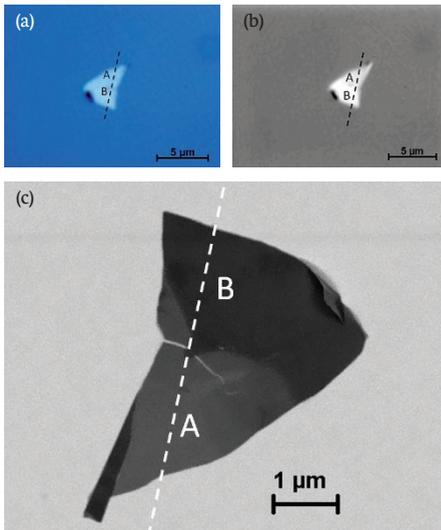

**Figure 5.** (a) Graphene layer stack observed in bright field and (b) with crossed circular polarization (red color channel). Optical phase profile measured along dashed line. (c) SEM of graphene layer stack on conductive silicon with 2nm native SiO<sub>2</sub> with In-lens Energy selective Backscatter-detector, extra high tension 1.5kV, grid bias 1.2kV, optical phase profile measure along dashed line

---

often sulphur. Scientists at the German microscopy company WITec have shown how a similar correlative microscopy workflow, combining light microscopy, Raman microscopy, SEM and AFM, can be used to study the TMDC tungsten disulphide ( $WS_2$ ). This involved using WITec's RISE microscope, which combines Raman microscopy and SEM in a single instrument.

## FURTHER INFORMATION

Bondia P, Jurado R, Casado S, *et al.* Hybrid nanoscopy of hybrid nanomaterials. *Small* 2017;13;1603784. (<http://dx.doi.org/10.1002/sml.201603784>)

Bradley RS, Withers PJ. Correlative multiscale tomography of biological materials. *MRS Bulletin* 2016;41(7):549–56. (<https://doi.org/10.1557/mrs.2016.137>)

Gelb J, Finegan DP, Brett DJL, *et al.* Multi-scale 3D investigations of a commercial 18650 Li-ion battery with correlative electron- and X-ray microscopy. *J Power Sources* 2017;357:77. (<http://dx.doi.org/10.1016/j.jpowsour.2017.04.102>)

Hauser M, Wojcik M, Kim D, *et al.* Correlative super-resolution microscopy: New dimensions and new opportunities. *Chem Rev* 2017. (<http://dx.doi.org/10.1021/acs.chemrev.6b00604>)

Merkle AP, Gelb J, Orchowski A, Fuchs J. X-ray microscopy: the cornerstone for correlative characterization methods in materials research and life science. *Microsc Microanal* 2014;20(Suppl 3):986. (<http://dx.doi.org/10.1017/S1431927614006655>)

Thomas C. Correlative microscopy of optical materials. *Imaging & Microscopy* Oct 13 2014. (<http://www.imaging-git.com/science/electron-and-ion-microscopy/correlative-microscopy-optical-material>)

Valdrè G. Correlative microscopy and probing in materials science. In Rickerby DG, Valdrè G, Valdrè U, eds. *Impact of electron scanning probe microscopy on materials research*. Dordrecht: Springer, 1999; 455–72. ([https://link.springer.com/chapter/10.1007/978-94-011-4451-3\\_22](https://link.springer.com/chapter/10.1007/978-94-011-4451-3_22))

ZEISS webinar: Multi-scale correlative microscopy of energy storage and conversion devices, June 2016. (<https://zeiss-microscopy.uberflip.com/h/i/242514217-multi-scale-correlative-microscopy-of-energy-storage-and-conversion-devices>)

